

Option Printer



A new optional printer is available for our laboratory autoclaves SX-300E/SX-500E/SX-700E, SX-300/SX-500/SX-700 to improve the overall quality of the product including safety control. It can print out the chamber temperature and pressure data, as well as operating conditions. The temperature data of sterilized items measured by an optional external temperature sensor can also be printed.

Supplementary information		
Product information	Error information	Information about changed settings
Model/Serial Number/Date and Time/Work; total number of operations.	When some error conditions occur during a run, information about the error is outputted.	When any setting has been changed during a run, the name of the changed item is outputted.

Data output image in each operating cycle

Time	Elapsed time since start of operation	TE	Temperature of sterilized items measured by an external temperature sensor
TC	Chamber temperature	P	Chamber pressure

```

==DATA=====
<Time> Time [hour:minute' second]
<TC> Chamber Temp. [°C]
<TE> External Temp. [°C]
<P> Chamber Press. [MPa]
- <Time> --- <TC> --- <TE> --- <P> ---
    
```

```

Temperature raising
00:00 19 0.000
00:03 20 0.000
00:06 23 0.000
:
    
```

```

Temperature & pressure raising
(00:22' 09)
00:24 105 0.025
00:27 113 0.065
    
```

```

Sterilizing & Count start
(00:29' 30 121 0.103)
00:30 121 0.103
00:31 121 0.105
    
```

```

Temperature & pressure reduction
(00:49' 34)
00:51 118 0.091
00:54 114 0.067
:
    
```

```

Temperature reduction
(01:06' 16)
01:09 98 0.000
    
```

```


Work completion.
(01:09' 17)
    
```

```

==DATA=====
<Time> Time [hour:minute' second]
<TC> Chamber Temp. [°C]
<TE> External Temp. [°C]
<P> Chamber Press. [MPa]
- <Time> --- <TC> --- <TE> --- <P> ---
Temperature raising
00:00 19 0.000
00:03 20 0.000
00:06 23 0.000
:
Temperature & pressure raising
(00:22' 09)
00:24 105 0.025
00:27 113 0.065
Sterilizing & Count start
(00:29' 30 121 0.103)
00:30 121 0.103
00:31 121 0.105
Temperature & pressure reduction
(00:49' 34)
00:51 118 0.091
00:54 114 0.067
:
Temperature reduction
(01:06' 16)
01:09 98 0.000
Work completion.
(01:09' 17)
=====
    
```

<Accessories included>

Thermal paper 1roll
Paper holder 1pc
Paper guide 1pc

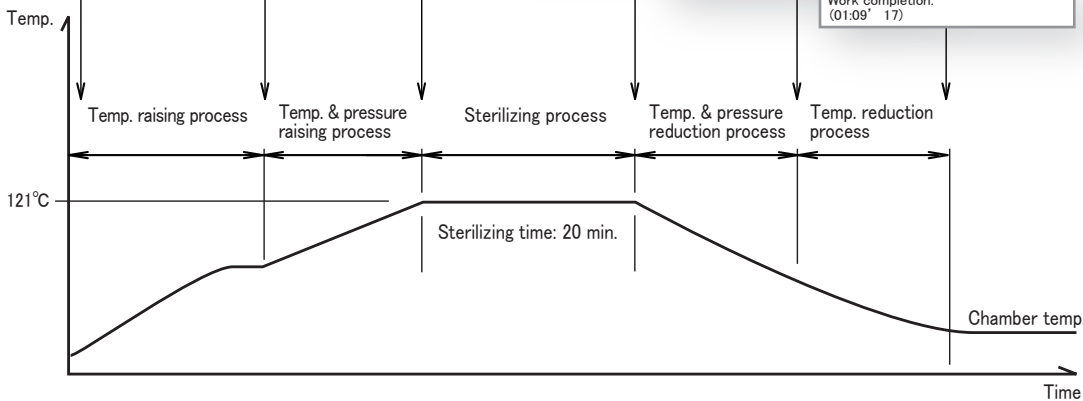

○ Specifications are subject to change according to product advancement.

Sales Office

TOMY DIGITAL BIOLOGY CO., LTD.

3-14-17 Tagara, Nerima-ku, Tokyo 179-0073, Japan
e-mail : info@digital-biology.co.jp
URL : http://www.digital-biology.co.jp
phone : +81-3-5971-8160
fax : +81-3-3970-6036

Manufacturer

TOMY KOGYO CO., LTD.

3-14-17 Tagara, Nerima-ku, Tokyo 179-0073, Japan